

## Easy and fast sequencing using safe barcode labels!



DNA barcode (orange) for

- Run24
- Run24Supreme

Request labels free of charge!



DNA barcode (green) for

- Run24Barcode

Please inquire for special discounts!



Primer barcode (yellow)

Request labels free of charge!

1.

Activate barcodes in myGATC

Welcome Sequencing Software ... and more Company myGATC

You are here: > myGATC > Barcode activation

**Activate barcodes**

In order to have the barcode numbers available in your myGATC online order form, you need to activate them first.

For barcode activation please enter batch (Batch Identification Number) below. These numbers are available on your labels (see the figure on the right side).

Batch number

BIN

**Activate**

2.


Choose activated barcode in your order form

b) define my sequencing jobs right here:

#	Template	Primer
1	Type: <input type="text" value="-"/> Name: <input type="text" value="new DNA name"/> DNA @ GATC Prepaid barcode: <input type="text" value="-"/> <ul style="list-style-type: none"> <li>TP0001</li> <li>MP0103</li> <li>JS1104</li> <li>JS1107</li> </ul>	Name: <input type="text" value="new Primer name"/> Primer @ GATC Sequence: <input type="text" value="Primer sequence"/> Barcode: <input type="text" value="-"/> Synthesis? <input type="checkbox"/>

No more jobs to define

3.

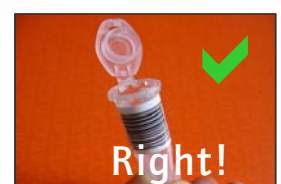

Please affix the label to the FRONT of the tube!